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**From its natural development to its processing, coal contains an amount of moisture which needs to be quantified.**

One method of quantifying moisture in coal is by measuring the weight loss of a sample after drying. The MFS is used for this drying process and has a small heated chamber to provide the lowest practical volume, ie, minimum free space, as required by the test standards.

A known mass of the coal is heated in a stream of nitrogen or air at a temperature between 105°C and 110°C and maintained at this temperature until its mass is constant. The moisture content is calculated from the loss in mass of the coal.

The ovens have a corrosion and oxidation resistant aluminium chamber which provides good temperature uniformity. The nitrogen or air flow can be adjusted by a flow meter fitted on the control panel and passes through a pre-heating chamber before entering the front of the work chamber.

The MFS/1 ISO operates with a regulated flow of moisturefree bottled nitrogen which removes the moisture released by the coal at 105°C. The MFS/1 ASTM operates with a regulated flow of air.

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