STOLP OVEN - V-L

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[Klik om video te bekijken](#)

Product Video: V-L soldering tube furnace

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- | Precise defined atmosphere with highest possible purity (6N or better)
- | Best possible vacuum
- | Designed for rapid heating and cooling
- | Certified safety management for flammable and toxic gases
- | Fully automatic operation
- | Movable quartz recipient
- | Data registratie voor kwaliteitsbeheer

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CIM, MIM, Quenching, drogen, gloeien, ontbinden, ontgassen, pyrolyse, rapid prototyping, sinteren, solderen, soldering, sublimering, synthese, temperen

STOLP OVEN - V-L

TECHNISCHE DETAILS

The soldering furnace is insulated with ceramic fibre with heating elements of CrFeAl wire. The top of the furnace hearth is closed and sealed with a ceramic fibre plug. The furnace has three heating zones controlled by the operating software to achieve the best possible temperature uniformity. The temperature of the soldering furnace is monitored and controlled by mantle thermocouples in each zone, in addition to, an over-temperature thermocouple.

The maximum temperature is limited to 1050 °C, which is the maximum possible temperature for the quartz tube under vacuum. The quartz tube is closed at the top end with the bottom end sealed to a flange at the loading plate. Metallic radiation shields are inserted to insulate the heat towards the loading plate. At the loading plate, the pumping unit is attached.

Located below the loading plate, a gas guiding tube is attached and positioned such that it is 300 mm above the ground plate to ensure gas flow is from the top to the bottom of the quartz tube. The gas outlet is attached to the vacuum flange. From this vacuum flange, several sample thermocouples can be positioned as required in the quartz tube. The furnace is operated by a fully, programmable operating system with a touch panel display for monitoring the process temperature and operating parameters.

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Opgelet

* at a height of 300 mm with a three zone furnace

www.carbolite.com/vl